

Single cell molecular characterization and PD-L1 expression analysis of model CTC using the RareCyte platform

Lance U'Ren, Yao Sun, Nolan Ericson, Tad George
RareCyte, Inc., Seattle, WA

RARECYTE

Background

Circulating tumor cells (CTC) are seeds for metastasis, the mechanism most responsible for cancer-related deaths. CTCs can be detected in peripheral blood, affording oncologists the attractive potential to phenotype a patient's cancer and to monitor changes in disease over time using a non-invasive procedure. Evaluation of CTCs are thus an attractive approach to understanding the potential response or non-response to anti-PD-1/PD-L1 therapies. CTCs are challenging to identify and characterize because they are exceedingly rare events within blood. RareCyte has developed the AccuCyte® - CyteFinder® system, which provides an exquisitely sensitive and repeatable workflow from blood collection to single CTC isolation. In a blinded fashion, we evaluated the sensitivity of RareCyte platform to detect model spike-in CTCs and correctly phenotype them based on PD-L1 expression and molecular characterization.

Sample preparation and analysis workflow

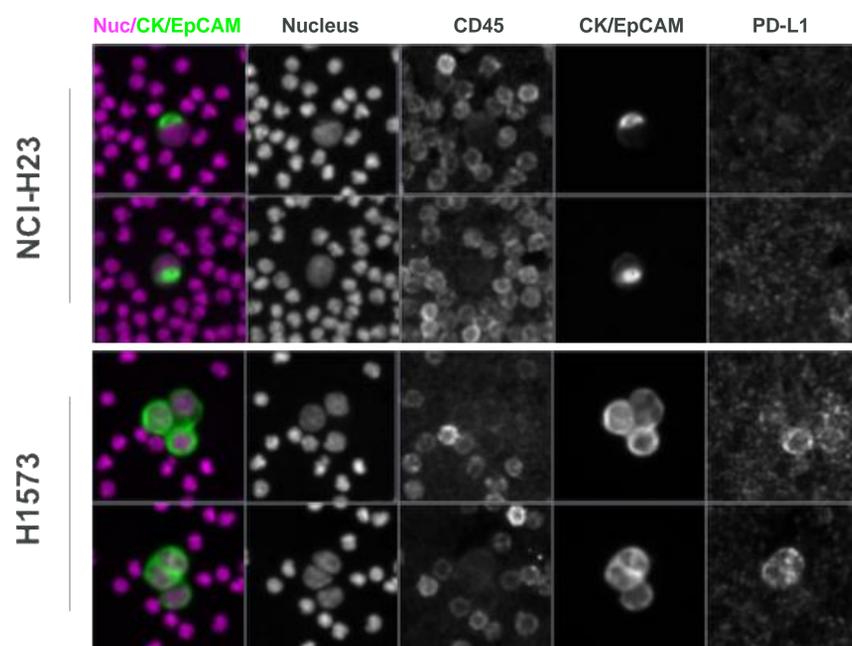
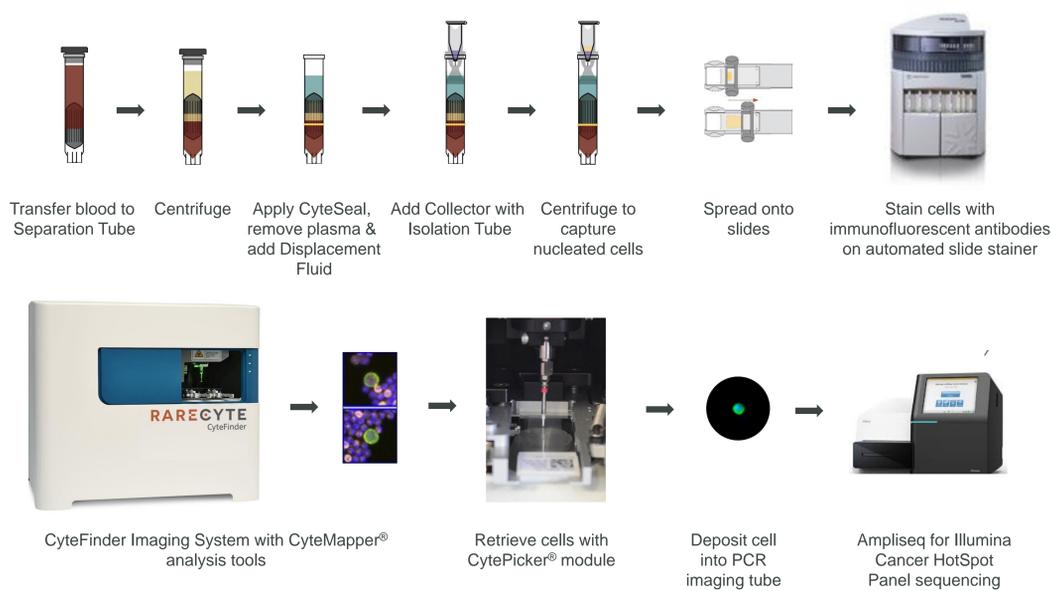


Figure 1: PD-L1 CTC spike-in models. 4-parameter analysis of cells detected using CyteFinder Imaging System

Sample	Cell Line	CTC Count [®] RareCyte	CTC Count [®] Pharma Partner	PDL1 MFI: Mean ± Std Dev	PDL1 MFI: Range	PDL1 MFI: Med (IQ Range)
1A	A549	81	106 - ND	154 ± 25	17 - 1597	84 (48 - 195)
2A	NCI-H358	86	118 - ND	364 ± 24	63 - 1327	322 (207 - 461)
3A	None	0	N/A	NA	NA	NA
4A*	NCI-H441	**4	23 - 5	898 ± 187	455 - 1362	887 (455 - 1260)
5A	NCI-H23	**16	26 - 11	148 ± 40	20 - 578	66 (53 - 227)
6A	MCF7	**7	21 - 16	60 ± 20	28 - 156	44 (31 - 83)
6B		**4				
7A	SK-MEL-28 [^]	0	16 - 9	NA	NA	NA
8A	H1573 ^{^^}	3	22 - 18	167 ± 53	129 - 228	144 (129 - 228)
8B		85				
9A	MDA-MB-231	7	12 - 9	312 ± 148	73 - 534	323 (202 - 400)
10A	NCI-H441	46	80 - ND	930 ± 405	67 - 1994	957 (661 - 1110)
11A	NCI-H23	7	9 - 9	92 ± 72	16 - 194	57 (34 - 154)
12A	SKBR3	14	17-10	887 ± 424	18 - 1636	979 (675 - 1098)

Table 1: PD-L1 CTC summary

® Cell count per 7.5 mL; RareCyte = CyteFinder; Pharma Partner = ViCell (pre-spike) - Calcein staining (post-spike)
* 4 of a total of 8 slides were stained/enumerated
** Clots visible around float for all samples processed on 180613; may affect CTC recovery
^ melanoma cell line that does not express EpCAM or CK and therefore not detected by the RareCyte panel used
^^ highly clumpy cell line

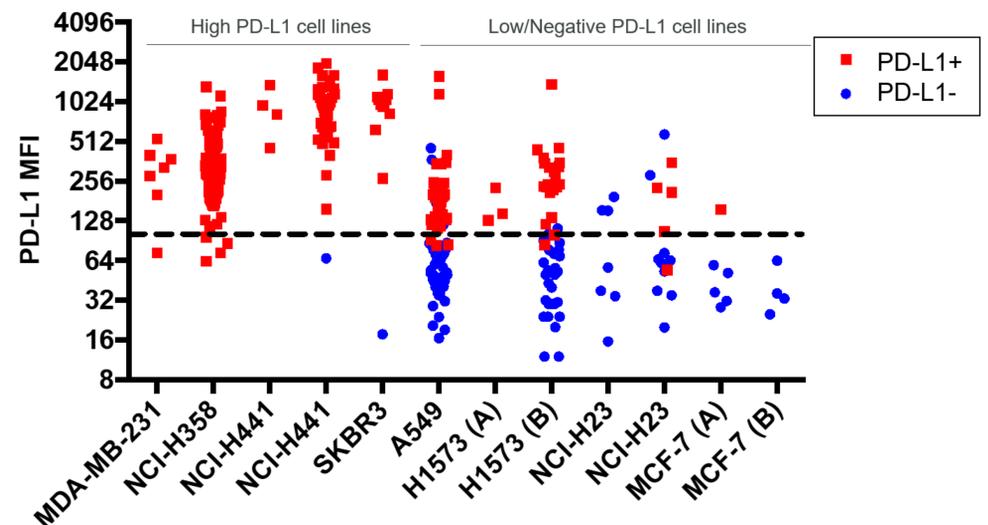


Figure 2: PD-L1 expression on CTC spike-in models. PD-L1 MFI (Mean Fluorescent Intensity) per pixel per cell using CyteFinder. CTC were visually stratified as PD-L1+ (■) or PD-L1- (●). Dashed line represents the 75th interquartile of all PD-L1- cells.

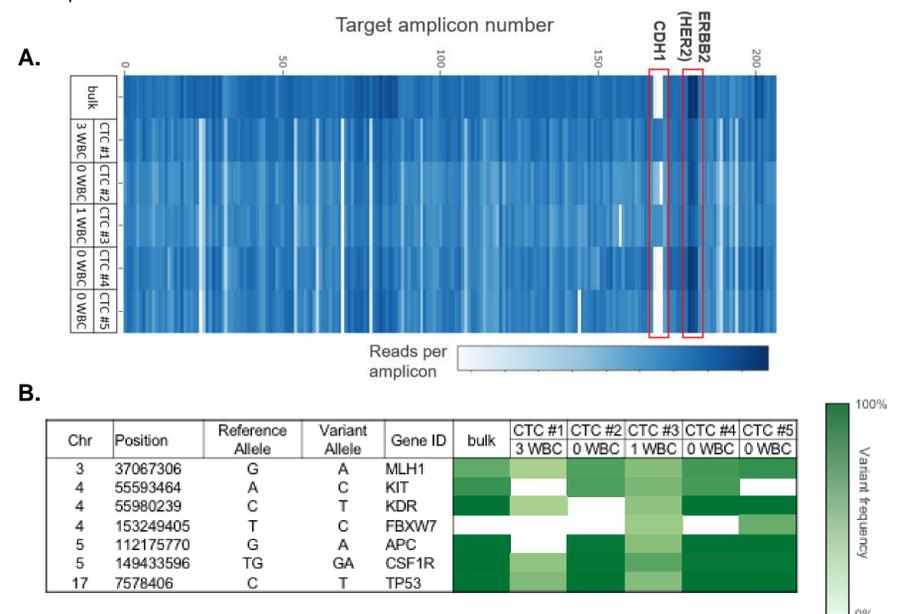


Figure 3: Targeted single cell DNA sequencing. A subset of spike-in cells were retrieved using the CytePicker Module and single-cell DNA sequencing was performed on a MiSeq (Illumina) using AmpliSeq™ Cancer Hotspot Panel (Illumina). A representative coverage heatmap displays the number of reads per target amplicon (A). Variant analysis was performed to identify cell line-specific mutations in individual cells (B).

Cell line	Expected mutation(s) [®]	Identified by single cell sequencing
A549	KRAS G12S STK11 Q37stop	✓ ✓
NCI-H358	TP53 deletion KRAS G12C	✓ ✓
MCF-7	PIK3CA E545K	✓
MDA-MB-231	TP53 R280K BRAF G464V* KRAS G13D*	✓
SKBR3	HER2 amplification CDH1 deletion MLH1 S406N KIT M541L TP53 R175H FBXW7 Y458C	✓ ✓ ✓ ✓ ✓ ✓

Table 2: Single cell targeted DNA NGS variant analysis summary.

® as determined by sequencing bulk genomic DNA isolated from cultured cells

* variants did not meet the detection threshold in single cells due to high numbers of co-retrieved white blood cells

Results

Upon sample unblinding, it was revealed that no CTCs were identified in no spike controls nor spikes of non-epithelial lines. Of the 8 epithelial lines tested, recovery frequencies of 7 were closely correlated with expected values, with the exception of the H1573 line, likely due to large cell clusters. PD-L1 expression correlated with expected results and mutations identified by sequencing matched published reports.

Conclusions

In a blinded study using 8 different cell lines as model CTC the RareCyte platform produced expected results for recoveries, PD-L1 phenotype and mutational analysis of retrieved CTCs. These results provide confidence in the utility of the RareCyte platform for clinical CTC enumeration, PD-L1 assessment, and single cell molecular characterization.